



Bedienungsanleitung • Mode d'emploi • Instruction

**KALTLICHTQUELLEN
SOURCES DE LUMIERE FROIDE
COLD LIGHT SOURCES**

MA 1100 / MA 1200 / MA 1300

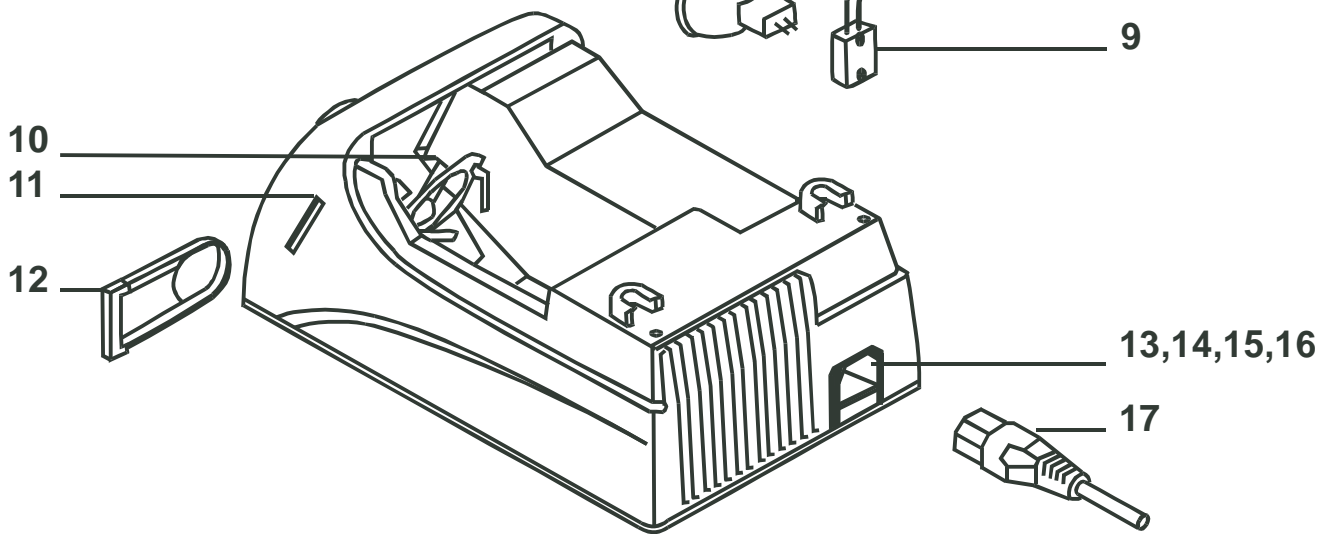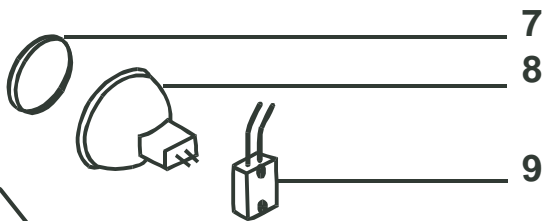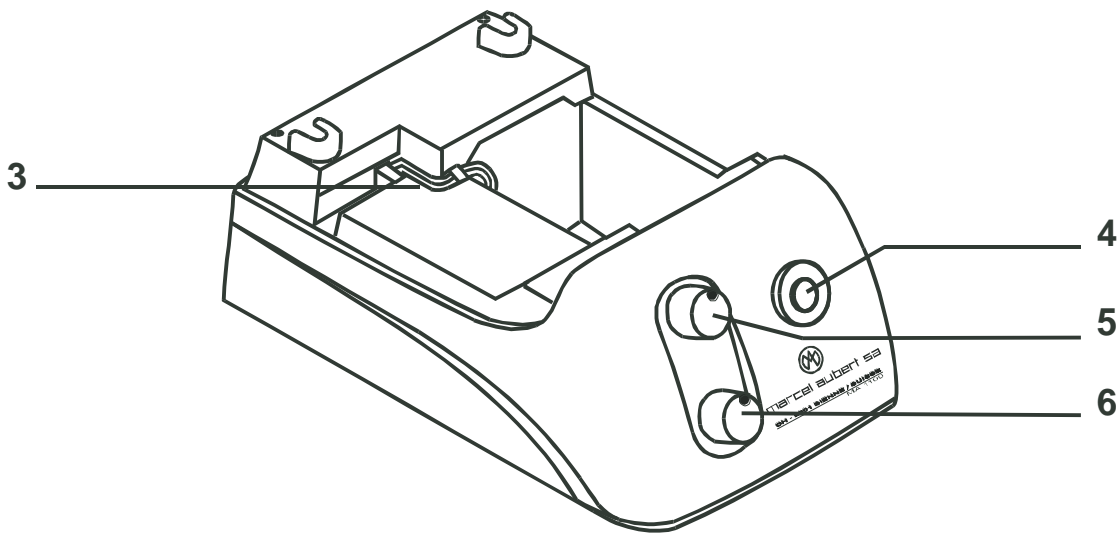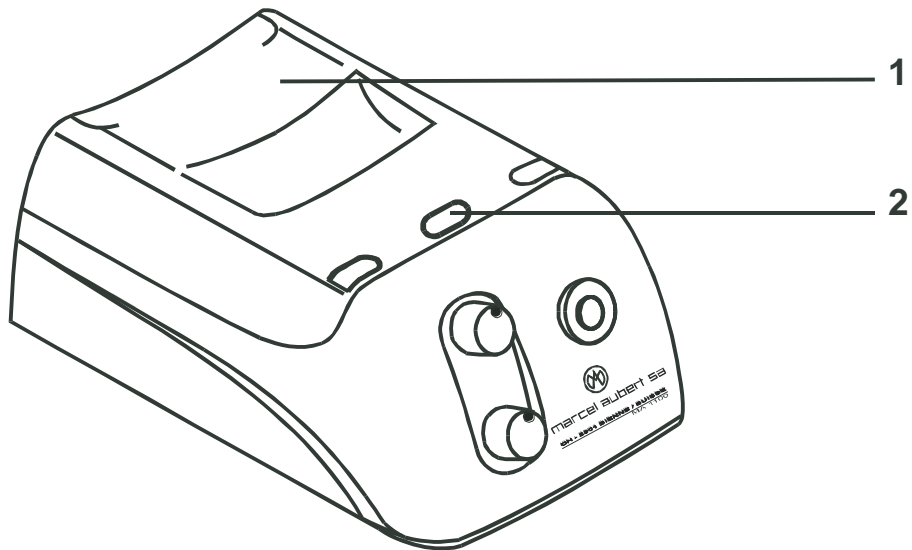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

Warning symbols

Caution: Risk of electrical shock!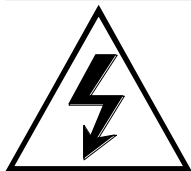

Caution: Dangerous area. Refer to accompanying documentation!

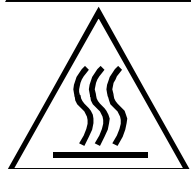Warning: Hot surfaces!

This section contains safety information, which must be observed strictly when using this device.

It is in your own interest to pay attention to all warnings on the device and in these operating instructions.

Target audience: Persons who work with this device or accessories of this device must read these operating instructions and the operator must ensure that the users understood the content of these instructions. This applies especially to the chapter entitled "Safety information".

Obligation to observe product: The user must immediately stop operation in case of changes of the operating behaviour or to safety-relevant components of the device (pull the power plug!) and inform the responsible person or the manufacturer.

Location of operating instructions: We recommend storing these operating instructions near the device to ensure quick access by operating staff.

Statutory provisions: National and local safety and accident prevention regulations and laws which are in force must be strictly observed in addition to the operational guidelines issued by the operating entity.

No technical modifications must be made to this device under any circumstances!

Caution! The lamp and parts in its immediate vicinity may be hot. The hot lamp and adjacent parts can cause burns upon contact. Before removing the cover and changing the lamp, the device must be allowed to cool off sufficiently (approx. 10-15 minutes). The power plug must be pulled first!

The halogen reflector lamp must be correctly placed for optimum function.

Never look directly into the lamp or the light guide when it is switched on; otherwise, you risk eye injuries!

Refer to specifications in the "Technical Data" section for binding operational limits.

Do not use this unit near water or in any space with excessive humidity. Do not place flammable materials on or near the unit at any time. Keep unit away from any sources of heat. The product has not been approved for operation in areas subject to explosion hazards! Not suitable for use with flammable mixtures (such as anaesthetics, oxygen or disinfectants....)!

Never block the ventilation openings of the instrument, otherwise it can overheat. Ensure that the free circulation of air is not impaired in any way.

Never spill liquid on the unit or insert objects into the unit! This could result in electric shock or damages to the unit.

The device may only be operated using the power supply voltage indicated.

Never open the device or any components unless instructed expressly to do so by these instructions.

Cleaning: Disconnect unit from the power supply before cleaning and only clean with a damp cloth. Never use combustible or flammable liquids. If fluids accidentally enter the device, unplug the power supply cable immediately and take the unit to the service location.

Maintenance: The operator must maintain all national laws and standards for regular inspections. An inspection must be performed not later than every 2 years by specialists who evaluate the safety of the device, recognize possible effects and dangers and can determine the conformity with guidelines and standards. These inspections should cover at least: Protective conductor inspection, leakage current inspection and visual inspection with cleaning and functional test

Spare parts: Only original spare parts may be used. If this is not done, it can lead to personal injury and material damage. Refer to the „Spare Parts“ section for component codes.

Liability: As the ultimate legal entity, the operating institution is responsible for ensuring the proper use of the device and for providing other operators with the necessary information and specifies the competencies required to operate the device.

Disposal: Observe all national environmental requirements and laws for the disposal of the device or its parts. Separate the components (e.g. by plastics, electronics, lamps, metals...) and dispose of them through national disposal companies!

Disconnect the power supply cable when the unit is not being used for an extended period of time. Use only original power cables and route the cables so that they cannot be pinched or separated. For safety reasons, only use a 3-wire power cable with protective conductor.

Repairs that are not described in this document must only be carried out by authorized workshops!

The manufacturer is not liable for any damage resulting from a failure to comply with the aforementioned instructions!

State of the art: This product constitutes state-of-the-art technology and employs recognized safety standards.

Select the future location based on the following criteria:

- No high humidity (see technical data).
- Normal room temperature (see technical data).
- Keep air vents free.
- Not suitable for use with flammable mixtures (such as anaesthetics, oxygen or disinfectants...).

Remove the device from the packaging and place it onto a sufficiently large and horizontal surface.

Do not yet connect the power plug.

Before initial operation, insert the halogen lamp. The device cover must be removed for this purpose.

- Press the release button (2).
- Remove the cover (1).

During the following operating steps, do not touch the quartz envelope of the inside of the reflector! This could lead to a premature failure of the lamp.

- Remove all transport packaging (e.g. carton for lamp socket).
- Fit lamp socket (9) on contact pins of the lamp without tilting.
- Insert lamp (8) into lamp holder (10) without tilting and push to the stop. If the lamp features an alignment feather, observe the correct engagement.
- Attach the cover (1) (self-engaging).
- Insert the fibre-optic light guide in the connection (4) to the stop. For fibre-optic light guides with feather pin, it must be inserted in the key to the stop. Ensure that the rotary knob (5) is in position "Off" (0-position).
- Insert the power plug.
- Activate the device by turning the rotary knob (5) clockwise.

The device may only be operated using the power supply voltage indicated. The specified power supply voltage can be found on the identification label at the bottom of the device. Only connect the unit to grounded sockets.

Keep air vents free!

Do not operate the device with open cover (1)!

The device has not been approved for operation in potentially explosive atmospheres!

Switch the device ON or OFF

- Operate the rotary knob (5).
- Position 0: Device is switched OFF.
- Position 1: Device is switched ON.

Adjust brightness (electrically)

- Turn the brightness knob (5).
- Turning the brightness knob clockwise continuously increases the lamp voltage and the brightness. Turning it counter clockwise reduces the brightness. Only adjust as much brightness as required because the useful life of the bulb decreases with increasing brightness.

Adjusting brightness with mechanical aperture

If you want to change the brightness without changing the color temperature, you should operate the mechanical hole matrix aperture. Turning the knob (6) clockwise increases the brightness up to the maximum. Turning it counter clockwise reduces the brightness up to approx. 1% of the maximum setting. These settings do not affect the life of the bulb. Of course, it is possible to operate the electrical brightness control, although the color temperature changes again. If both brightness controls are used, it results in a control range of 100% to approx. 0.1%.

Replacing the slide-in filter

- Remove the slide-in filter mount (12) from the slide-in filter guide (11).
- Slide the desired filter with mount into the slide-in filter guide until it engages.

Color and fluorescence filters can be used here for specific applications.

Use only original filters; otherwise, a thermal destruction of the filter and the device cannot be ruled. Note that the filter and the entire holder may become very hot during operation. Let the device cool down sufficiently before removing the filter!

TROUBLESHOOTING

- First, check whether the power cable is inserted at the light source and in a working power socket.
- Ensure the light source is switched on.

No light

Cause	Fault elimination
Defective bulb:	Replace bulb
Defective fuse:	Replace primary circuit fuse.
Thermal protector has responded:	Keep air vents clear, instrument restarts after a short time.
Defective lamp socket:	Replace lamp socket
Defective switched-mode power supply:	Return device for repair

Insufficient light

Cause	Fault removal
Bulb remains at the minimum setting (dimmed) despite moving the electrical brightness knob:	Return device for repair
Fan failure:	Return device for repair

Note: A defective bulb can also cause a fan standstill or a speed reduction.

REPLACING DEFECTIVE PARTS

Use only original spare parts! If this is not done, it can lead to personal injury and material damage. Repairs that are not described in this document must only be carried out by authorized workshops!

Before opening the device, always pull the power plug!

Do not touch the quartz envelope or the inside of the lamp reflector! This drastically shortens the life of the bulb. Grasp the bulb only at the outside of the reflector or close to the contact pins. Always let hot bulbs and their environment cool off sufficiently.

Replacing the bulb

Light source and bulb must be cooled off sufficiently before changing the bulb!

- Switch off the device (5) and pull the power plug (17).
- Let the defective bulb cool off (min. 10 minutes).
- Press the release button (2).
- Remove the cover (1).

During the following operating steps, do not touch the quartz envelope and the inside of the reflector!

- Pull up the defective halogen lamp (8) together with the lamp socket (9) from the lamp holder (10).
- Pull the defective bulb (8) out of the lamp socket (9).
- Insert the new bulb in the lamp socket (9) without tilting.
- Insert lamp (8) into lamp holder (10) without tilting and push to the stop. If the lamp features an alignment feather, observe the correct engagement.
- Attach the cover (1) (self-engaging).

Replacing the lamp socket

As soon as the lamp socket (2) shows burn-off signs between the contacts, it must be replaced. As a guiding value, a replacement of the lamp socket is recommended after 5 replaced bulbs.

- The procedure is identical to that described in the chapter "Replacing the bulb" up to the item "Remove the defective bulb (8) from the lamp socket (9)".
- Release 2 screws in the clamping bead (16) and pull out the old lamp socket.
- Pull the new lamp socket through the loop, insert it in the clamping bead and tighten the screws.
- The additional procedure is identical to that described in the chapter "Replacing the bulb" starting with the item "Insert new bulb in the lamp socket (9) without tilting".

Replacing the heat-absorbing filter (MA 1200 and MA 1300)

Gas the heat-absorbing filter only at the outer diameter! Do not contaminate the surfaces. Light source and heat-absorbing filter must be cooled off sufficiently before replacement. The heat-absorbing filter continues to be very hot after being switched off!

- Switch off the device (5) and pull the power plug (17).
- Press the release button (2) and remove the cover (1).
- Completely remove the broken heat-absorbing filter (7).
- Insert the new heat-absorbing filter in the spring guide
- Attach the cover (1) (self-engaging).

Replacing the primary fuse

Use only original fuses according to the chapter "Technical data".

- Switch off the device (5) and pull the power plug (17).
- Pull out the fuse tray (14) (snap lock).
- Remove defective fuses (15) and replace them with spare fuses (16).

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