
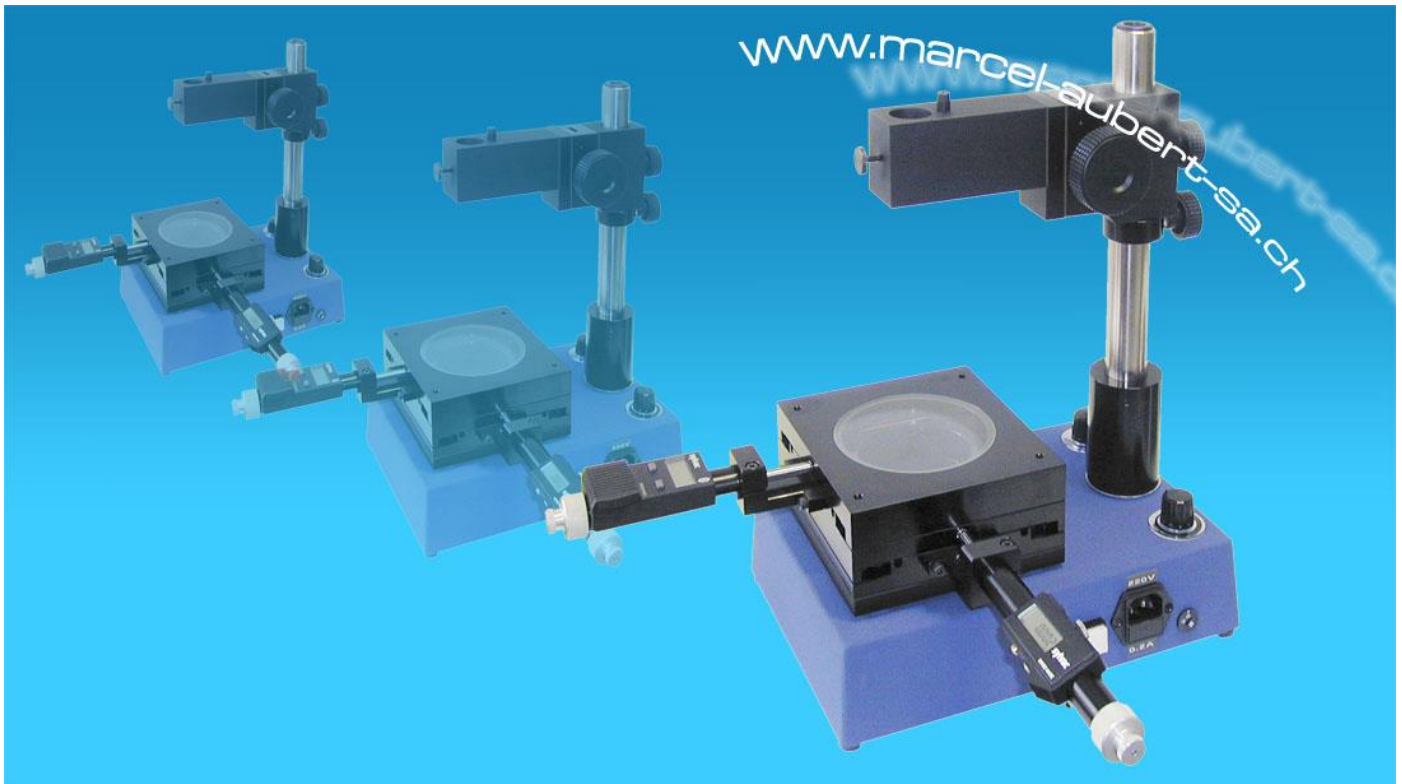
**Bedienungsanleitung • Mode d'emploi • Instruction**

**STATIV • STATIF • STAND**

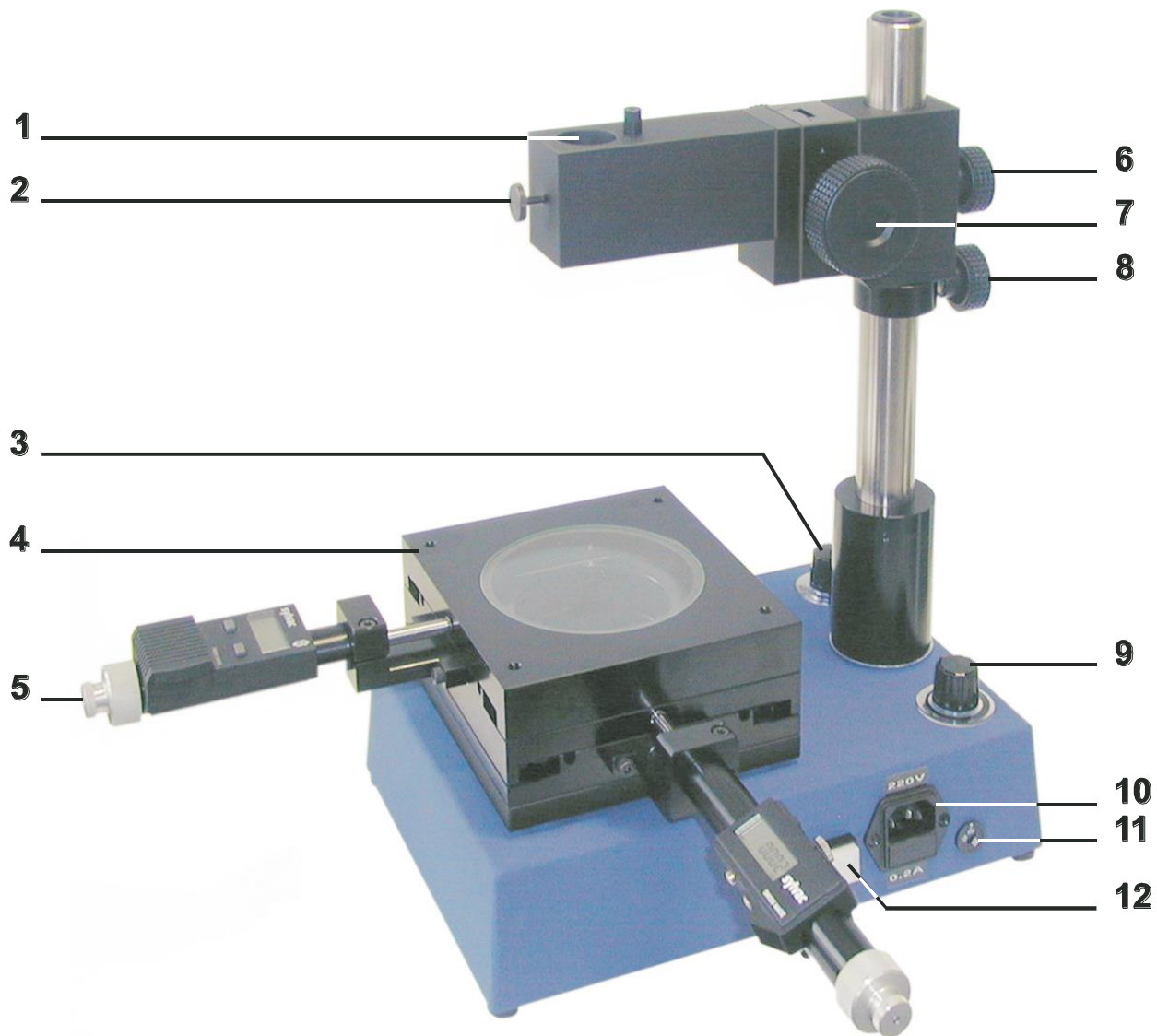**MA 331ES**

---

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---

1	Optic holder
2	Blocking screw (tube)
3	Intensity regulator of the transmitted light illuminator
4	Co-ordinate table
5	Micrometer heads
6	Blocking screw for the manual vertical adjustment
7	Vertical adjustment
8	Stop collar for vertical adjustment with blocking screw
9	Intensity regulator for incident light illuminator
10	Power supply line
11	Voltage switch 110/220V
12	Main switch

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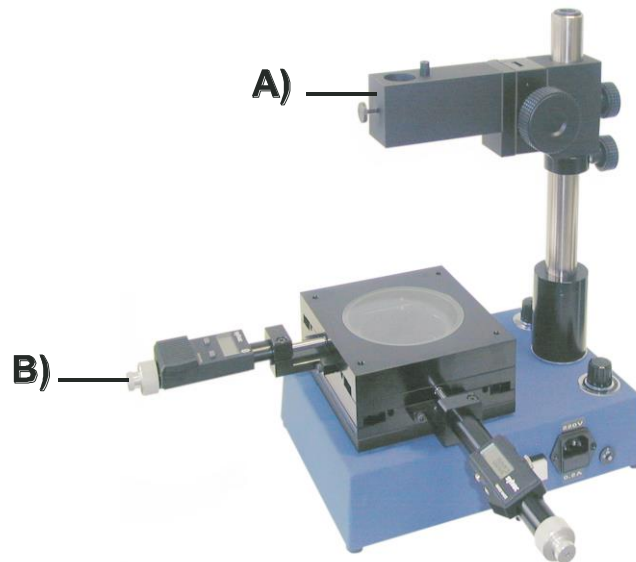
### **General security measures and notes:**

- Ensure that power supply corresponds to the unit voltage.
- Ensure that the power supply cord is in proper condition before using the unit.
- Take out the mains power supply before attempting to carry out any work.
- Do not touch the socket, bulb or lamp while the unit is functioning (heat).
- Some of the unit components, such as the table glass, screen, lens and others, can have sharp edges and should be operated with due care.
- Do not look at the light during long periods.
- Please pay the necessary attention by every displacement of the axis.
- Repairs should only be carried out by authorized or qualified staff.

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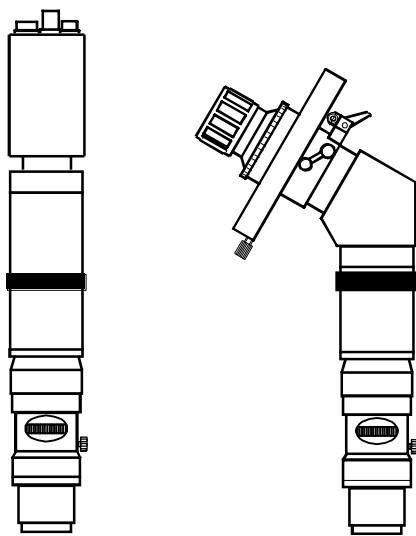
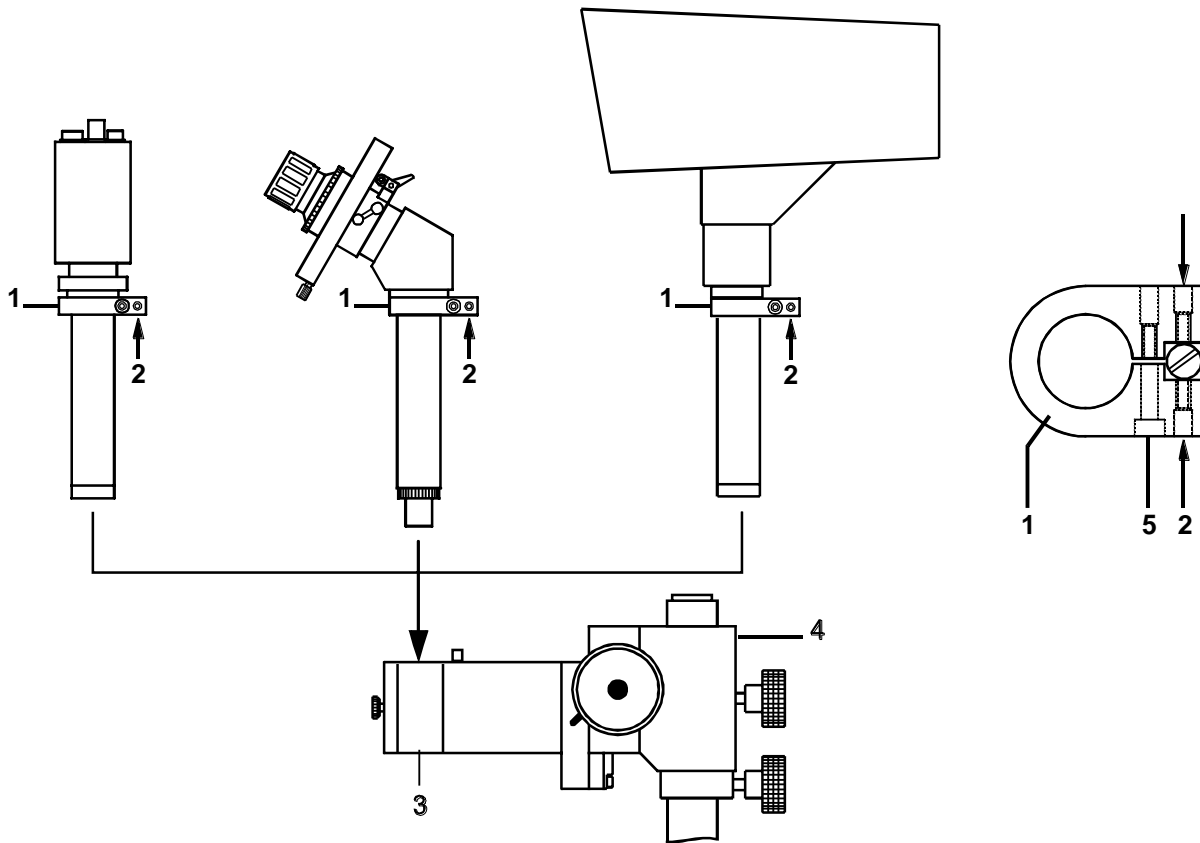
## Security measures for the transport

**Please ensure that**

- A) ... the unit is not lifted up by the vertical holder of the base.
- B) ... the unit is not lifted up by the micrometer heads.  
... generally, that all retaining screws are fixed correctly.

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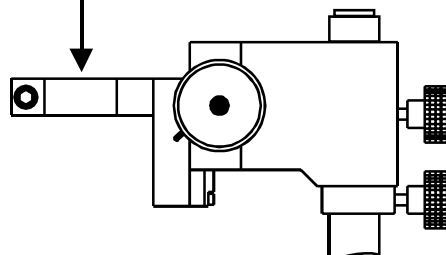
**The zoom-optic:**

See the instruction "Mounting of the zoom-optic"

1	<b>Verdrehsicherungsbride</b>	<b>Bride de positionnement</b>	<b>Fixing clamp</b>
2	<b>Schraube (rechts)</b>	<b>Vis (à droite)</b>	<b>Screw (right side)</b>
3	<b>Optikaufnahme</b>	<b>Support - optique</b>	<b>Optic holder</b>
4	<b>Vertikalträger des Grundgerätes</b>	<b>Support vertical de la base</b>	<b>Vertical holder of the base</b>
5	<b>Blockierschraube</b>	<b>Vis de blocage</b>	<b>Blocking screw</b>

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The optic (video, microscope or projector) has been carefully centred and checked before dispatch. It was removed for simplify the packing.

### **Mounting of the optic:**

- Mount the optic with fixing clamp (1) in the holder (3).
- Draw the lateral screw (2) which is positioned on the right side at the back of the fixing clamp. The optic has to come back at its correct position.
- For the correction of eventual alignment faults, adjust with the screw (2) and the screw which is positioned opposite.

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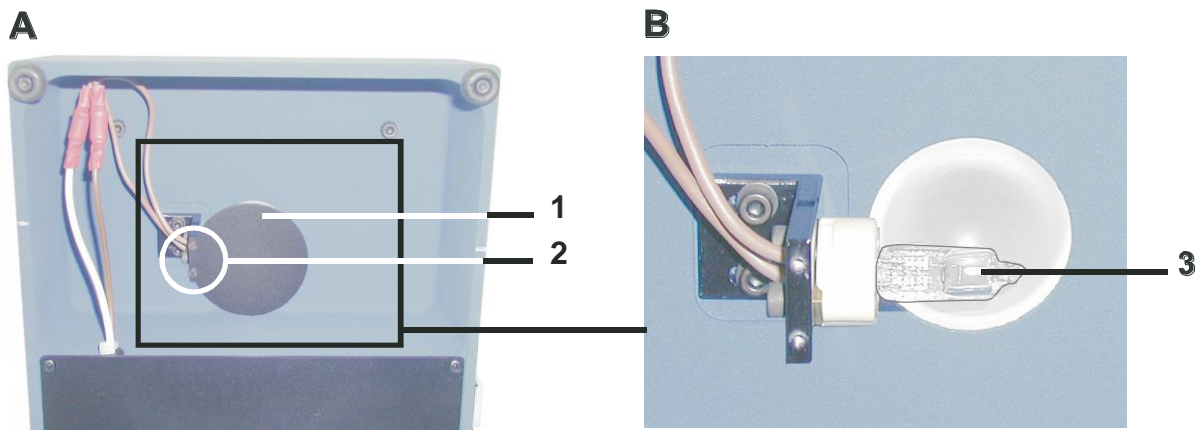
**Only valid for units with halogen lighting**

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## Renewal of the halogen-bulb

**Attention: Before manipulate the halogen-bulb let it cool down !**

- Disconnect plug.
- Unscrew the lateral screw which is positioned on the right side at the back of the fixing clamp (see "Mounting of the optic").
- Take down the optic (video, microscope or projector) from the base.
- Tip over backward the base, see picture A.
- Screw off the 2 screws (2) and take down the dazzle protection (1).
- Set a new halogen-bulb in (3), see picture B. **(Never grip the halogen bulb with bare fingers).**
- Mount the dazzle protection (1) and screw on the 2 screws (2).
- Set up the base and mount the optic (see "Mounting of the optic").
- Connect plug.
- Regulate the brightness with the intensity regulator for transmitted illuminator.  
**(Attention: Danger of dazzling!)**

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