
ZUBEHOER  
ACCESSORY  
ACCESSOIRE

Bedienungsanleitung • Mode d'emploi • Instruction  
**Titan View**

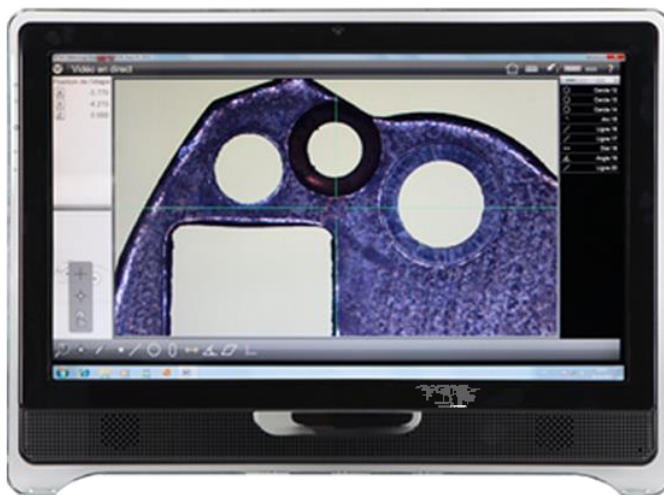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
## Before Power up

### Electrical connection

Line voltage 100 V~ to 240 V~ (-15 % to +10 %)  
Line frequency 43 Hz to 63 Hz

---

### Power connector wiring

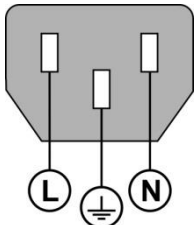L : Line voltage (brown)

N : Neutral (blue)

 Earth ground (yellow/green)

---

### Danger of electrical shock!

- Do not open the enclosure.
  - Never use 3-wire to 2-wire adapters.
  - The insulating jacket of the cable shouldn't have any defects
- 

### Caution

General safety precautions must be followed when operating this electrical equipment. Do not observe these precautions could damage the equipment or injure personnel. Safety rules vary within companies. If a conflict exists between the content of this guide and the rules of a company using this system, the more stringent rules should be applied.

---

## Functions

- 1) Light control ( Intensity setting / Mode selection)
- 2) Mode control
- 3) Optional :Push-button (Light sector selection)
- 4) Master switch

## General purpose

This light source series was developed to perform a quick and high precise setting of the light intensity.

**Option:** Sector illumination with the ring light.

- Turn the device on by pushing the master switch (4) (illuminated indicator= device is working).
- Set the luminous intensity with the potentiometer (1).
- Push the potentiometer (1) to select the setting mode:
  - **Coarse** mode ( control (2) = green) to quickly reach the interesting light range,
  - **Fine** mode (control (2) = orange), approximately 10 times finer, to precisely adjust the image contrast.

The coarse setting is active when the device starts.


The last luminous intensity used is automatically saved before the device is switch down (a short stabilisation time <1 sec is required).

---

## Specifications for ring light with sectors

Some ring lights can illuminate samples in a more directional way (especially useful for surface observation) and can be controlled by sectors.

Use the push-button (3) to select the most adapted light solution. Eight configurations are available.

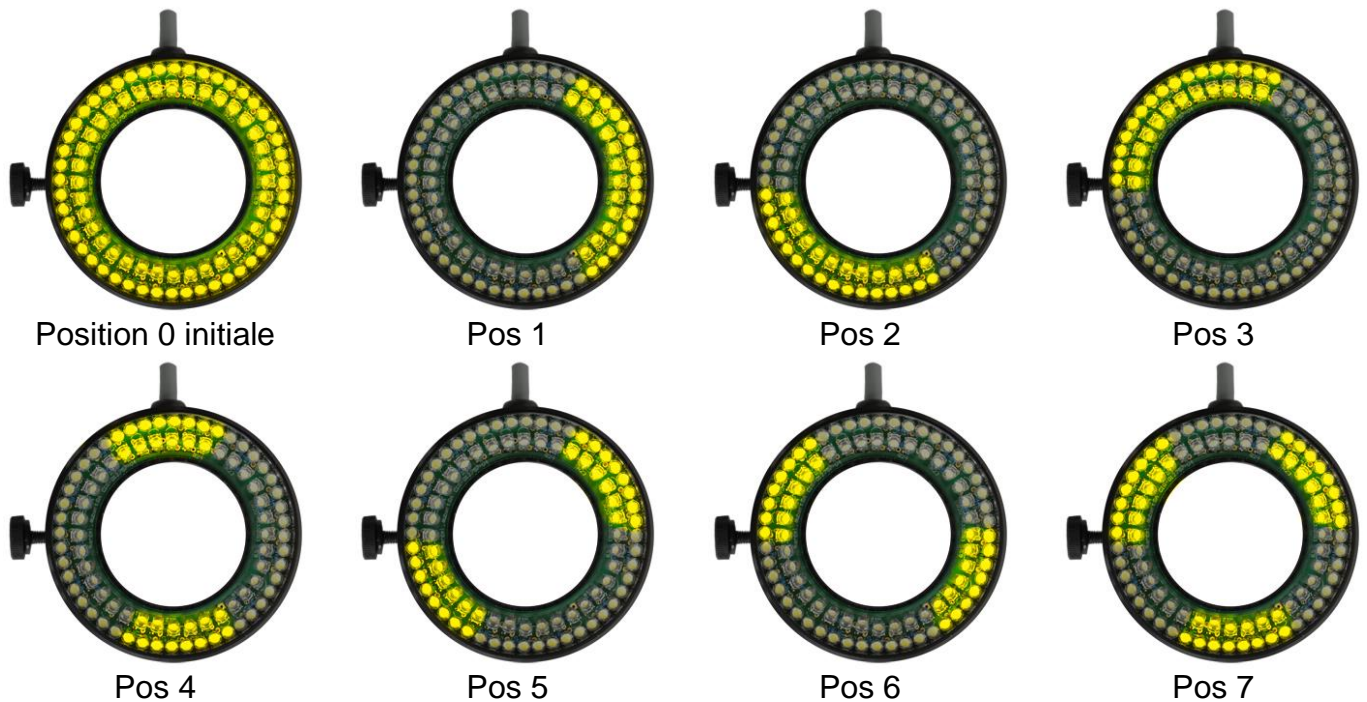The coarse setting and the default **position 0** are active when the device starts.

The last light intensity used on each sector is also automatically saved for each configuration.

## Troubleshooting

Selection of sectors unresponsive	Turn off the power, wait minimum 3 seconds, and turn on the power. (Master switch No. 4).
Selection of sectors responds slowly	Check that the potentiometer is not positioned between two notches. (Light control No. 1).
Potentiometer does not change the intensity.	Faulty potentiometer (factory return).
Potentiometer does not change in coarse or fin mode	Turn off the power, wait minimum 3 seconds, and turn on the power. (Master switch No. 4).
On / off or Master switch does not light up.	Defective switch (factory return), or switch can be replaced by a qualified technician (Art. No. BPL-017712)

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