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
Important information before powering on

Electrical connection

Line voltage 100 V~ to 240 V~ (-15 % to +10 %)

Line frequency 43 Hz to 63 Hz

Power connector wiring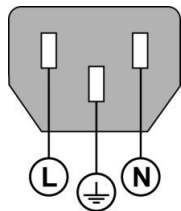

L : Line voltage (brown)

N : Neutral (blue)

 Earth ground (yellow/green)

Danger of electrical shock!

Caution


- Do not open the enclosure.
- Never use 3-wire to 2-wire adapters.
- The insulating jacket of the cable shouldn't have any defects

General safety precautions must be followed when operating this electrical equipment. Do not observe these precautions could damage the equipment or injure personnel.


Safety rules vary within companies. If a conflict exists between the content of this guide and the rules of a company using this system, the more stringent rules should be applied.

Use of incident light sectors


Our 6-sector ring light incident illuminators (Ø25 or Ø40) offer the possibility to illuminate samples selectively by controlling light by sectors. This makes it possible to highlight different contrasts according to the angle or the surface state of the part to be observed. To change the sector position, press button (**3**) for MA 356-041 and **5**) for MA 356-042) successively until you obtain one of the 8 available configurations. The incident light sectors are reset to 0 (all sectors switched on) during ignition so as not to disturb users.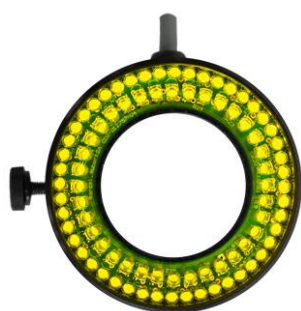


Pos 0 « initial »

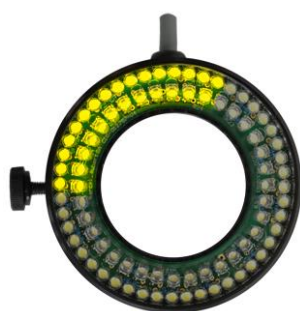
Pos 1

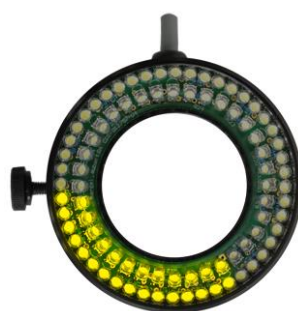
Pos 2

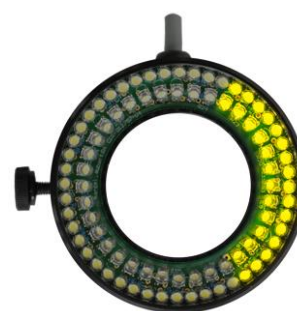
Pos 3

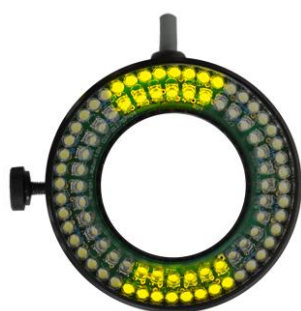
Pos 4

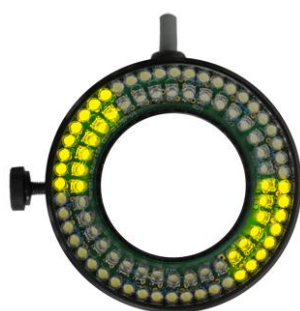Pos 5

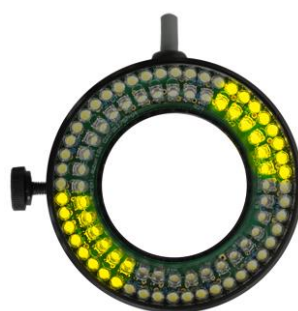Pos 6

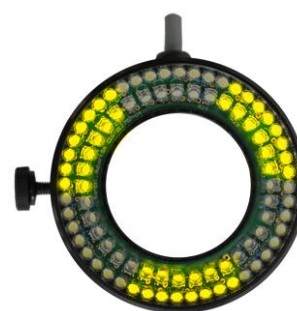Pos 7

MA 356-041-...

Descriptions

- 1) Light control (Intensity setting / Mode selection)
- 2) Mode control
- 3) Optional :Push-button (Light sector selection)
- 4) Master switch

General purpose

This light source series was developed to perform a quick and high precise setting of the light intensity.

Option: Sector illumination with the ring light.

- Turn the device on by pushing the master switch (4) (illuminated indicator= device is working).
- Set the luminous intensity with the potentiometer (1).
- Push the potentiometer (1) to select the setting mode:
 - **Coarse** mode (control (2) = green) to quickly reach the interesting light range,
 - **Fine** mode (control (2) = orange), approximately 10 times finer, to precisely adjust the image contrast.


The coarse setting is active when the device starts.

The last luminous intensity used is automatically saved before the device is switch down (a short stabilisation time <1 sec is required).

Troubleshooting


Selection of sectors unresponsive	Turn off the power, wait minimum 3 seconds, and turn on the power. (Master switch No. 4).
Selection of sectors responds slowly	Check that the potentiometer is not positioned between two notches. (Light control No. 1).
Potentiometer does not change the intensity.	Faulty potentiometer (factory return).
Potentiometer does not change in coarse or fin mode	Turn off the power, wait minimum 3 seconds, and turn on the power. (Master switch No. 4).
On / off or Master switch does not light up.	Defective switch (factory return), or switch can be replaced by a qualified technician (Art. No. BPL-017712)

Descriptions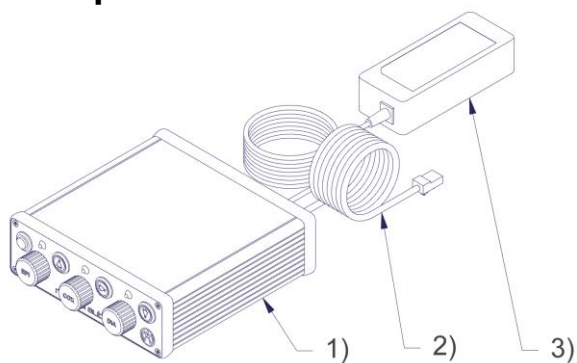


- 1) Illuminators control box
- 2) Ethernet cable
- 3) Power supply 100/240V
- 4) Indicator light of the illuminator mode in use.
- 5) Incident illuminator sector selector switch (circular lighting from the top)
- 6) Incident illuminator intensity regulator
- 7) Incident illuminator On/Off switch
- 8) Coaxial illuminator intensity regulator (lighting passing through optical axis)
- 9) Coaxial illuminator On/Off switch
- 10) Transmitted illuminator sector selector switch (lighting from the bottom)
- 11) Transmitted illuminator On/Off switch
- 12) Ethernet connector to plug cable 2) into computer (only for software that can control illuminators)
- 13) « Power jack » connector for power supply 3)
- 14) Transmitted illuminator connector
- 15) Coaxial illuminator connector
- 16) Incident illuminator connector

MA 356-042-01 « Opensys »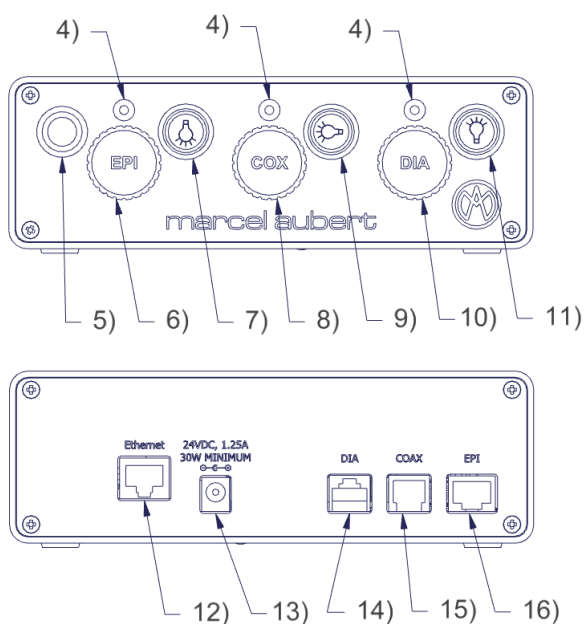

MA 356-042-02 « Rotatif »

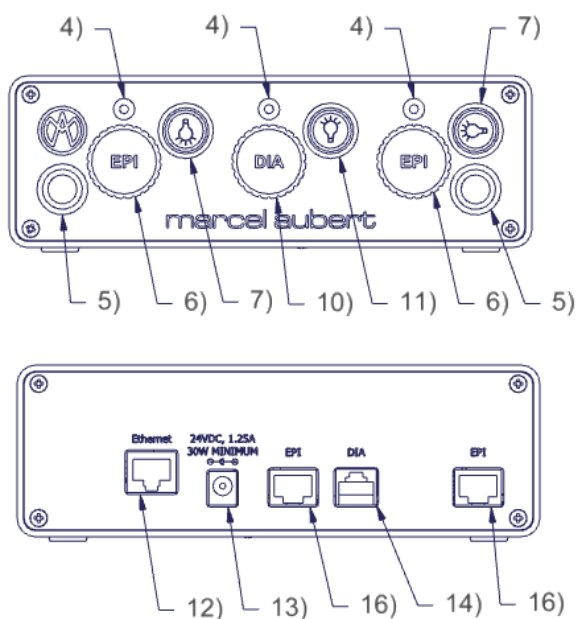General use

- The unit is always on. It was designed to allow a precise and fast intensity adjustment of the different LED illuminators.
- The box automatically saves the last lighting intensity used by the operator (a short delay <1 second is required before the device can be powered off).
- You can independently turn the illuminators on or off. To do this, push the On/Off switch of the corresponding illuminator ; **7**), **9**) or **11**).
- The light intensities can be adjusted by turning the knobs **6**), **8**) and **10**)
- Each lighting has a coarse and fine mode. To switch from one to another, press the corresponding intensity button, ie; **6**), **8**) or **10**).
- Coarse mode (indicator light **4**) = green), to quickly reach the desired illuminating range.
- Fine mode (indicator light **4**) = orange), about 10x plus more sensitive, for optimum contrast control of the image.
- By default, when the device is switched on, the coarse mode is active.

Troubleshooting

Sector selector switch 5) no longer responds.

Disconnect the power supply **13**), wait for at least 3 seconds and switch on the power again.

All lights are off on the housing **1**).

Power supply **3**) is defective. You can order this article : ALM-021462.

Buttons **6**), **8**) or **10**) do not change the intensity of the illuminator.

The knob is defective, the box **1**) must be returned to the factory.

Buttons **6**), **8**) or **10**) do not change between coarse / fine mode.

Disconnect the power supply **13**), wait for at least 3 seconds and switch on the power again.

On/Off switches **7**), **9**) or **11**) do not work anymore.

The switch is defective, the box **1**) must be returned to the factory.

For any other case, please contact us.

Tel: +41 (0)32 365 51 31
Fax: +41 (0)32 365 76 20
Email: info@marcel-aubert-sa.ch

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